

## Instructions for: HFN6

### *HF Acid Eater Neutralizer & Absorbent*

Read and follow all instructions and warnings on the product packaging  
**AVOID ALL CONTACT WITH ACIDS EXCEPT AS DIRECTED.**

#### **Assessment**

Identify the material spilled: Assess the toxicity, flammability, or other hazardous properties of the material, the size, type, amount and location of the spill.

#### **Preparation**

Identify and obtain your personal protective equipment. This might include (but is not limited to) safety goggles, facemask, dust mask, gloves, apron, booties or more.

#### **Clean Up**

1. Stop the source of the spill. Notify surrounding personnel of the spill, isolate and limit access to the immediate area where the spill has occurred.

2. Gently apply the HF Acid Eater Absorber & Neutralizer around the edges of the spill, building a dike or barrier to contain it.

3. Then, slowly pour the absorber, working from the outside of the spill inward, until the liquid is completely covered.

4. Allow HF Acid Eater™ Absorber & Neutralizer a few minutes to completely neutralize and absorb the spill. The product will form into a solid mass making it easy to clean up.

5. Prepare the disposal bags. Gently sweep up the material and place in the disposal bags.

6. Clean and rinse personal protective equipment with water before re-using.

7. Dispose of all used articles according to local, state and federal regulations.

